

# Primed T-Cell Activation: Antigen-Presenting Cell Co-Culture Method

---

## FC17

This method provides a general procedure for reactivating antigen-experienced (primed or memory) T-cells prior to staining for intracellular cytokines, activation markers, or cell proliferation. T cells are co-cultured with an antigen-presenting cell, in this case, dendritic cells that have been pulsed with an antigen of interest.

This protocol provides a general staining procedure for use with Bio-Rad reagents. Specific recommendations are provided on product datasheets, and those instructions should always be used in conjunction with the product- and batch-specific information provided with each antibody vial. A certain level of technical skill is required for the successful design and implementation of these techniques; these are guidelines only and may need to be adjusted for particular cell types or applications.

### Reagents:

- Cell culture medium
- Dendritic cells from an appropriate source, such as murine bone marrow cells cultured in 10 mg/mL GM-CSF
- Antigen of interest, such as a synthetic peptide
- Lipopolysaccharide (LPS)
- Required when measuring proliferation:
  - CytoTrack Cell Proliferation Assay Kit (catalog #1351202, 1351203, or 1351205) or CFDA-SE Cell Proliferation Assay Kit (#1351201)
- Phosphate buffered saline, 10x (PBS) (#BUF036A)
- Required when staining cell surface antigens:
  - Staining Buffer (#BUF073)



## Method:

1. Pulse dendritic cells with antigen (50–200 µg/mL) by incubating overnight in the presence of LPS (100 ng/mL). Antigen should be titrated to obtain reproducible results.  
**Note:** Non-pulsed dendritic cells should be included as a negative control.
2. Wash the dendritic cells twice in cell culture media. Resuspend cells in cell culture media and adjust to  $1 \times 10^6$  cells/mL.
3. Isolate T cells and resuspend in media at  $1 \times 10^6$  cells/mL.
4. **Optional:** For proliferation studies using CytoTrack Cell Proliferation Assay or CFDA-SE Cell Proliferation Assay, incubate cells with the dye following the recommended protocol, or see protocol **FC18**, Measuring Cell Proliferation Using CytoTrack, a Cell-Permeable Dye.
5. Co-culture the dendritic cells and T cells at increasing ratios, for example, 1:1, 1:5, and 1:10 (DC:T cells) to obtain a range of stimulation.
6. Incubate cells in a humidified 37°C, 5% CO<sub>2</sub> incubator for 6–96 hr as required.
7. Centrifuge at 300–400 x g for 5 min and discard supernatant.
8. Perform staining of cell surface antigens (protocol **FC4**, Direct Immunofluorescence Staining of Surface Epitopes of Cells and Whole Blood) or intracellular staining (protocol **FC7**, Direct Immunofluorescence Staining of Intracellular Antigens: The Leucoperm Method) as required along with associated wash steps.  
**Note:** Cytokine staining requires protein transport inhibitors such as brefeldin A and monensin. See protocol **FC9**, Direct Immunofluorescence Staining of Intracellular Cytokines in Whole Blood, for more details.
9. Acquire samples on a flow cytometer.

## Notes:

- Other types of antigen-presenting cells may be used but could require pulsing with different amounts of antigen and different incubation times
- All antibodies should be titrated prior to use to ensure that the optimal concentration is used
- Appropriate controls should be included; for example, an unstimulated control and incubation with non-pulsed dendritic cells

Visit [bio-rad-antibodies.com/flow](https://www.bio-rad-antibodies.com/flow) for more flow cytometry resources.

Bio-Rad is a trademark of Bio-Rad Laboratories, Inc. All trademarks used herein are the property of their respective owner.  
© 2026 Bio-Rad Laboratories.



**Bio-Rad**  
Laboratories, Inc.

Life Science  
Group

Website [bio-rad.com](https://www.bio-rad.com) USA 1 800 424 6723 Australia 61 2 9914 2800 Austria 00 800 00 24 67 23 Belgium 00 800 00 24 67 23 Brazil 55 11 3065 7550 Canada 1 800 361 1808 China 86 21 6169 8500 Czech Republic 00 800 00 24 67 23 Denmark 00 800 00 24 67 23 Finland 00 800 00 24 67 23 France 00 800 00 24 67 23 Germany 00 800 00 24 67 23 Greece 30 210 7774396 Hong Kong 852 2789 3300 Hungary 00 800 00 24 67 23 India 91 124 4029300 Israel 000 800 00 24 67 23 Italy 00 800 00 24 67 23 Japan 81 3 6361 7000 Korea 82 080 007 7373 Luxembourg 00 800 00 24 67 23 Mexico 52 55 5488 7670 The Netherlands 00 800 00 24 67 23 New Zealand 64 9 415 2280 Norway 00 800 00 24 67 23 Poland 00 800 00 24 67 23 Portugal 00 800 00 24 67 23 Russian Federation 7 495 721 14 04 Singapore 65 6415-3170 South Africa 27 21 531 7504 Spain 00 800 00 24 67 23 Sweden 00 800 00 24 67 23 Switzerland 00 800 00 24 67 23 Taiwan 886 2 2578 7189 Thailand 662 651 8311 United Arab Emirates 971 4 818 7300 United Kingdom 00 800 00 24 67 23